



Human Induced Pluripotent Stem Cells (from Normal, Male, Peripheral Blood Mononuclear Cells (PBMC))

Product Information

Catalog Number	ASE-9215
Description	Applied StemCell, Inc. provides Control Human Induced Pluripotent Stem (iPS) cells at low passages (p15). These pluripotent cells were generated from normal peripheral blood mononuclear cells (PBMC) using a Sendai virus method. This method allows the expression of human transcription factors (<i>OCT4</i> , <i>SOX-2</i> , <i>KLF4</i> , and <i>c-MYC</i>) that initiate the reprogramming process. The resulting human iPS cells (hiPSCs) were selected using morphological criteria without the use of either fluorescent markers or drug selection. These iPS cells have been tested for expression of pluripotency markers (Oct-4, Sox-2, Nanog, and SSEA-4; Figure 2) and for normal male karyotype (Figure 3). The ASE-9215 control human iPSC line can be used for CRISPR/Cas9 genome editing and differentiation to somatic lineages <i>in vitro</i> . Detailed protocols for thawing, culturing under feeder-free conditions, and cryopreservation of these iPS cells are provided.
Reprogramming Method	Sendai Virus
Passage #	p15
Tissue	Peripheral Blood Mononuclear Cells (PBMC)
Age	21 (Average)
Sex	Male
Race	Caucasian
Clinical information	Normal
Quantity	1x10 ⁶ cells/vial
Shipping	Dry ice
Storage and Stability	Store in liquid nitrogen freezer immediately upon receipt. This product is stable for at least 6 months from the date of receiving when stored as directed.
Quality Control	Each lot of human iPS cells has been tested for growth and viability following recovery from cryopreservation, morphology, immunohistochemistry for pluripotency markers (Oct-4, Sox-2, Nanog, and SSEA-4); karyotyping; and for the absence of bacteria, fungi, mycoplasma (CoA available upon request).
Safety Precaution	PLEASE READ BEFORE HANDLING ANY FROZEN VIALS. Please wear appropriate Personal Protection Equipment (lab coat, thermal gloves, safety goggles)

Applied StemCell, Inc.

521 Cottonwood Dr. #111, Milpitas, CA 95035

Phone: 866-497-4180 (US Toll Free); 408-773-8007 Fax: 408-773-8238

info@appliedstemcell.com www.appliedstemcell.com

Copyright 2022, Applied StemCell, Inc. All rights reserved. This information is subject to change without notice.

and a face shield) when handling frozen vials. Please be aware that the following scenario can occur: Liquid nitrogen can leak into the vials when the vials are submerged in liquid nitrogen. Upon thawing, the liquid nitrogen returns to the gas phase, resulting in a dangerous build-up of pressure within the vial. This can result in the vial exploding and expelling not only the vial contents but also the vial cap and plastic fragments of the vial.

Warranty and Liability

Applied StemCell (ASC) does not extend any expressed or implied warranties. ASC does not provide any warranties of merchantability or suitability for specific product use. Upon purchase, ASC does not provide any warranty that the use of the product will not infringe any patent, copyright, trademark, or other proprietary rights. The customer acknowledges that the product may be part of a patent application or covered by patent rights in one or more countries. No licenses, expressed or implied, are provided when product is purchased. No license or rights to use a third-party patent, technology, or intellectual property will be included in the purchase of the ASC product. The customer is responsible for obtaining the necessary permissions based on their intend use of the product from third parties that may have a proprietary interest in the product.

Liability Statement for Customers Receiving the ASC Product: Customer agrees that they will be held responsible for any damages, claims, costs, and expenses resulting from any injury (including death), damage, or loss that may be caused by the use or distribution of the ASC product purchased or any derivatives to the extent permitted by law.

Indemnification: Customer agrees to hold harmless Applied StemCell from any damages, claims, costs, or expenses that may result from use of the product and/or derivatives, including injury (including death), damage, or loss.

Restricted Use

This product is for research use only and not intended for human or animal diagnostic or therapeutic uses.

Characterization of iPSC Line ASE-9215

Morphology Image

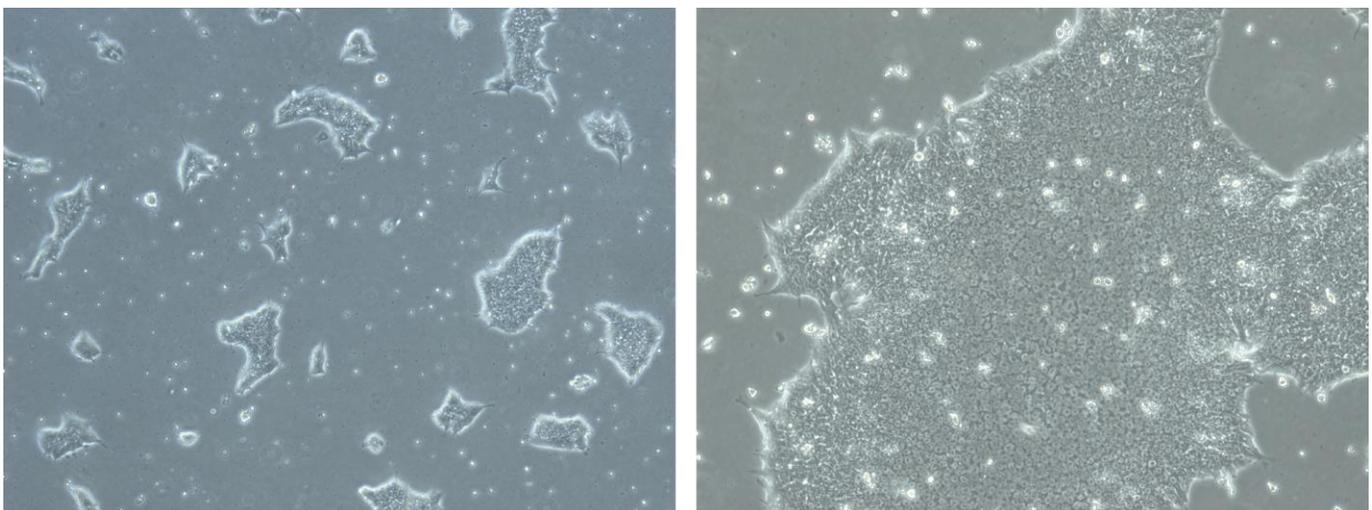


Figure 1: Morphology images of the ASE-9215 iPSC line captured 24 hours and 72 hours post thaw. The control human iPSCs were generated from normal Peripheral Blood Mononuclear Cells (PBMC) using a Sendai virus method. Feeder-free reprogramming media was used. The image on the left was taken 24 hours after thawing and the image on the right was captured 72 hours after thawing.

Pluripotency Marker Analysis

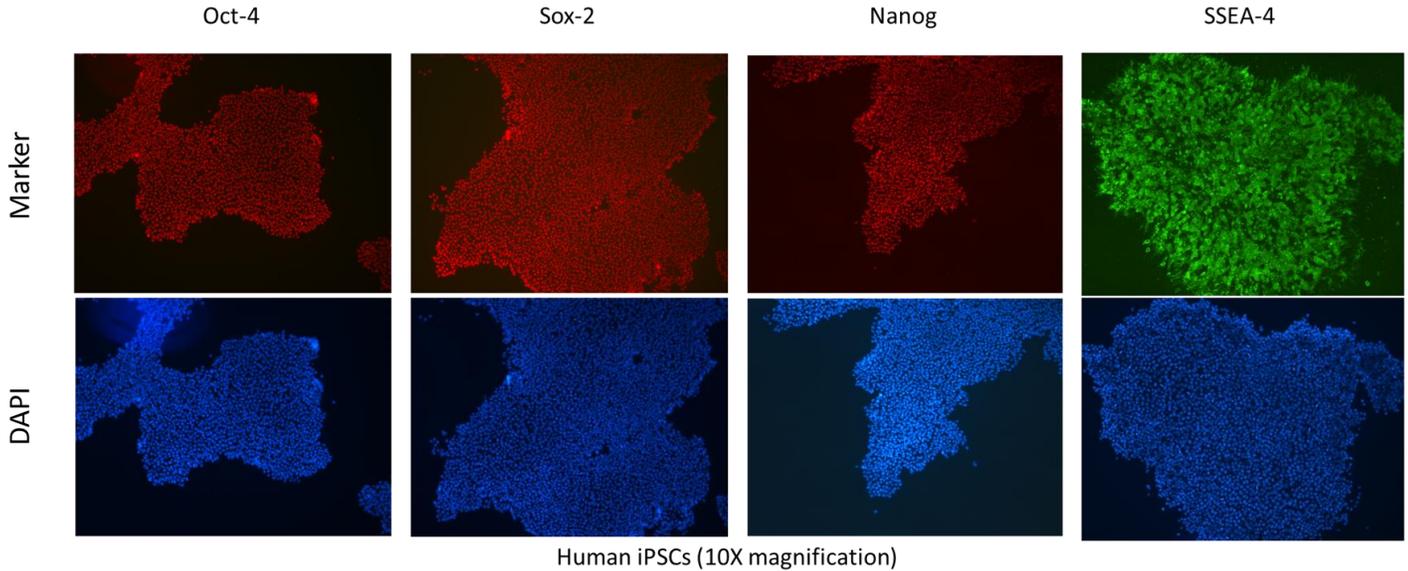


Figure 2: Expression of pluripotency markers. The ASE-9215 iPSC cell line expresses common iPSC biomarkers (top row: Oct-4, Sox-2, Nanog, and SSEA-4). Bottom row: Corresponding DAPI nuclear staining. All images were taken at 10X magnification.

Karyotype Analysis



Figure 3: Karyotype analysis to rule out genetic aberrations. Cytogenetic analysis was performed on twenty G-banded metaphase cells from human cell line ASE-9215 p5. Seventeen cells demonstrated an apparently normal male karyotype, and three cells demonstrated non-clonal chromosome aberrations, which are most likely artifacts of culture.

Media and Material Required but not Provided

- mTeSR-Plus Medium, StemCell Technologies, Cat# 05825
- Matrigel®, Corning, Cat# 354277
- ROCK Inhibitor (Y-27632), Sigma-Aldrich, Cat# SCM075
- CryoStor® CS10, StemCell Technologies, Cat# 7930
- 0.5M EDTA in PBS, ThermoFisher Scientific, Cat# 15575-020
- PBS, Life Technologies, Cat# 14190136
- 100 µg/mL Primocin™, Invivogen, Cat# ant-pm-2 or 1% Penicillin- Streptomycin, Life Technologies, Cat# 15140
- Cell Scraper, VWR International, Cat# 75799-938
- Corning® CoolCell® FTS30, Corning®, Cat# 432006

Protocol

Feeder-free culture conditions

1. Thawing human iPSCs using a feeder-free protocol

- 1.1 Prepare Matrigel® coated 6-well plates in advance, following vendor's instructions.
- 1.2 Prepare 6 mL of mTeSR-plus medium + 10 µM Rock Inhibitor.
- 1.3 Add 1 mL of mTeSR-plus medium + Rock Inhibitor in one well of a Matrigel®-coated plate. Prepare 2 wells for each vial of frozen cells.
- 1.4 Bring the cryovial on dry ice to the tissue culture room.
- 1.5 Quickly thaw the iPSCs in a 37°C water bath, by gently shaking the cryovial continuously until only a small piece of ice remains.
- 1.6 Wipe the cryovial with a paper towel sprayed with 70% ethanol and place it into a biosafety cabinet.
- 1.7 Add 9 mL of mTeSR-plus medium to a 15 mL conical tube.
- 1.8 Using a 1 mL pipette transfer the cells to the 15 mL conical tube dropwise while swirling the conical tube.
- 1.9 Centrifuge the cells for 3 min at 200 RCF and at room temperature.
- 1.10 Aspirate off the medium, and add 4 mL mTeSR-plus medium + Rock Inhibitor.
- 1.11 Gently flick the conical tube to resuspend the cells and transfer them to the 2 wells of the Matrigel® plate using a 5 mL serological pipette.
Note: Prepare Matrigel plates the day before and not more than 3 days prior to thawing. Before transferring the cells, the Matrigel must be aspirated.
- 1.12 Place the plate in the incubator and move the plate back-and-forth and side-to-side, twice to spread the clumps evenly in the wells.
- 1.13 Medium can be changed 2 days after thawing. After the 2 days, if the number of attached colonies is still low (less than 5% confluency), change half of the medium (aspirate 1 mL and add 1 mL of fresh medium).
- 1.14 Once colonies are stabilized, change the medium daily. Usually, within 1 week the cells are ready to be split.
- 1.15 When the ASE-9215 hiPSC colonies are big or close enough to merge, the cells need splitting/ passaging.

2. Passaging/ splitting human iPSCs using EDTA

- 2.1 Aspirate the medium from the hiPSC culture.
- 2.2 Wash once with 1 mL of 0.5 mM EDTA (in PBS).
- 2.3 Aspirate the EDTA and add 1 mL of 0.5 mM EDTA (in PBS) per well of a 6-well plate and incubate the cells for 3 min in a 37°C incubator.
- 2.4 Observe the cells under a microscope until the cells at the edge of the colonies start to separate and round up.
- 2.5 With the cells still attached, aspirate the EDTA and add 1 mL of mTeSR-plus medium.
- 2.6 Scrape the cells from the bottom of the well until the colonies are floating; pipette up and down 2-3 times (with the 1000 µL pipette set to 800 µL) to break the colonies in small clumps.
Note: Pipetting up and down three times is enough to break the colonies into clumps of optimal size.
- 2.7 Transfer the desired dilution to the wells of the new Matrigel®-coated plate (usually around a 1:3 splitting ratio).
Note: Prepare Matrigel plates (or any other basement matrix) not more than a week prior to passaging cells.

- 2.8 Place the plate in the incubator and move the plate back and forth and side to side, twice to spread the clumps evenly in all the wells.

Note: hiPSCs are passaged as clumps of 50-200 cells, rather than single cells. Cells need to be passaged before the colonies are large enough to merge with one another.

3. Cryopreserving human iPS cells

- 3.1 Label the cryovial as needed, based on 1 vial per well of a 6-well plate, and pre-chill them in a 4°C freezer.
- 3.2 Aspirate the medium from the hiPSC culture.
- 3.3 Wash once with 1 mL of PBS.
- 3.4 Aspirate the PBS and add 1 mL of 0.5 mM EDTA (in PBS) per well of a 6-well plate and incubate the cells for 3-5 minutes in a 37°C incubator.
- 3.5 Observe the cells under a microscope. After 1-2 minutes the cells at the edge of the colonies will start to separate and round up.
- 3.6 Once round colonies are observed, aspirate the EDTA (even if it hasn't been 5 minutes of incubation).
- 3.7 Add 1 mL of cold freezing medium, CS10 and scrape the cells from the bottom of the well until the colonies are floating.
Note: The freezing medium must remain at 4°C until usage.
- 3.8 Pipette the cells once or twice to break up any big clumps before transferring the suspension to a cryovial.
- 3.9 Transfer the cell suspension into the pre-chilled and labeled cryovial.
- 3.10 Place the cryovial in a CoolCell® Freezing Container or in a Styrofoam rack at - 80°C overnight, and transfer to liquid nitrogen the next day.